Tel: 400-968-6088 Fax: 010-56371281/82

Http://www.solarbio.com

中性树胶使用说明书

Neutral balsam

货号: G8590

规格: 100ml

保存:室温干燥保存,有效期2年

产品说明:

显微技术上使用的优良封藏剂,与加拿大树胶性质相似,用途相同,效果较好。中性树胶与加拿大树胶相同,是一种天然树脂,经提炼与中和后溶解在透明剂中成60%溶液,折光率与玻璃近似,干燥后凝结成透明无色的固体,没有收缩变黄、纹裂的弊病,没有像加拿大树胶有年久变黄切片氧化退色等缺点。切片标本有苏木精、曙红、番红、固绿、洋红染色的,保藏数年颜色无变化。许多整装标本如吸血虫、涤虫等寄生虫,用中性树胶封固,均能取得优良效果。

目前中性树胶广泛使用在生物学,医学作为生物切片封藏剂,以及光学琉璃仪器的粘固剂。

使用说明:

封片方法很简单,在载玻片中央滴 2-3 滴树胶液 (注意要滴在一起否则容易有气泡),然后把盖玻片慢慢放上去,轻轻挤压,树胶液会向外侧扩散至整个镜片,通风晾几个小时后用棉球粘上透明剂或者松节油把镜片边缘多余的树胶擦洗干净。然后再放在通风处晾干,封片可常温保存。

注意事项:

透明剂有毒,避免与皮肤或眼睛接触。封片时注意通风。

相关产品:

P1010 0.01M PBS 干粉, PH7.2-7.4

C1010 0.01M 柠檬酸钠缓冲液 pH6.0

X1020 胰蛋白酶消化液 0.1%

T1080 20×*TBS*

DA1010 DAB 显色试剂盒

PR1100 BCIP/NBT 显色试剂盒

P2100 10×多聚赖氨酸

S2100 抗荧光衰减封片剂

相关文献:

- [1] Xinrui Zhao,Fufu Zheng,Yizhuo Li,et al. BPTF promotes hepatocellular carcinoma growth by modulating hTERT signaling and cancer stem cell traits. Redox Biology. July 2018. (IF 7.126)
- [2] Yuzhe Sun, Jiafeng Li, Zhendong Wei, et al. Proteomic and bioinformatic analysis of condyloma acuminata: mild hyperthermia treatment reveals compromised HPV infectivity of keratinocytes via regulation of metabolism, differentiation and anti-viral responses. International Journal of Hyperthermia. March 2019. (IF 3.589)

注: 更多使用本产品的文献请参考索莱宝官网。